

Multipurpose Electron Microscope Specimen Box Instructions



1. Record card is stored underneath the box

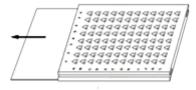


2. When moving lid and plastic cover pieces only four holes are exposed at once

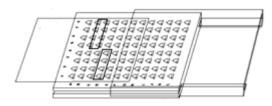



3. The tweezers insertion slot grips the grid edge



4. Slot grids should best be gripped by curved tweezers. Mesh grids can be gripped by straight or curved tweezers.

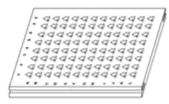

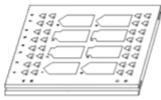

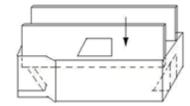

- 5. MEM-96grid contain 96 grids (left) MEM-32grid-8block contain 32 grids and 8 blocks (middle).
- 6. Small semi-thin section can be stored by sliding into the block cavity (right).