## **Technical Data Sheet**

## **Diamond Knife Wash Instructions**

## #<u>70605</u>

- 1. Put the knife wash on a Magnetic Stirrer. (make sure there is NO heat)
- 2. Fit the Diamond Knife into the holder and tighten the screw making sure that the screw meets the rear of the boat.
- 3. Fill the wash to the indicated line with distilled or deionized water (you may add two drops of very mild liquid detergent).
- 4. Put the lid on the container (be careful when doing).
- 5. Turn the magnetic stirrer on at a low speed for 5 minutes.
- 6. Remove lid and take knife out.
- 7. If desired and/or you used detergent, rinse the knife with distilled water.
- 8. Either let the knife dry naturally or if you prefer you can dry it with a foam rod (polystyrene) dipped in absolute ethanol as usual.





