New Packaging syringe Lead Citrate EM stain with Reynolds Formulation

Adapted

TEM Stainer RMC QG3100 Automate grids TEM staining


Reference :

ref # 11300-QG



The syringe 30 ml AirLess is used to dispense the lead citrate solution modified after Reynolds.

30 ml Lead citrate with formulation Reynolds

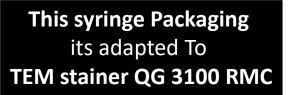

-6

7

9

_ml/cc -12

30 ml LEAD CITRATE STAIN Reynolds <mark># 11300-QG</mark> for TEM Stainer RMC QG3100

2. Install the lead citrate syringe in the syringe holder by removing the brass block from the holder, and inserting the syringe with its outlet downwards.

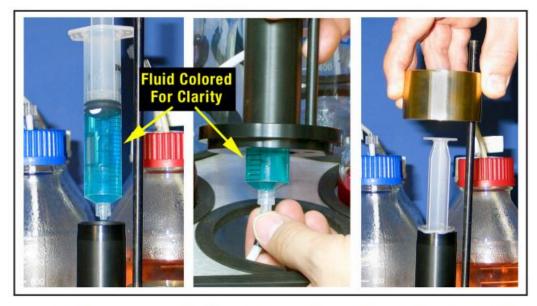


Figure 2: 1) Insert syringe in holder. 2) Connect tubing. 3) Gently lower brass block.

3. Connect the outlet of the syringe to the Stain 2 tube.

The connection with the tube must be done before the brass block is installed again.

30 ml LEAD CITRATE STAIN Reynolds ref # 11300-QG





4. After connecting the tube, gently guide the black rod through the hole in the brass block until it rests on the plunger.



Figure 3: Syringe and holder ready for use.







